

1. Installation of USB2EPP cable drivers for Windows

a) Attach USB2EPP cable to spectrometer. **b)** Attach UP5V power adapter to spectrometer. Green LED illuminates. **c)** Attach USB2EPP cable to available USB port on PC. Insert StellarNet CDROM and allow WinXP to search for a suitable driver.

If not WinXP then specify USB2EPP cable driver location on CDROM.

For Win00/XP = "\SWDrivers-USB2EPP\Win2k"

For Win98/Me = "\SWDrivers-USB2EPP\Win9x"

Use Device Manager to verify "StellarNet USB2EPP cable" is listed under USBDEV device. On WinXP, right click on MyComputer then select Manage, then select Device Manager.

If "...cable - start" appears, then disconnect cable from PC and remove UP5V then repeat **b)** and **c)** above, to re-install cable driver and verify proper driver installation.

Note: If you have the time, it is smart to repeat the USB2EPP cable installation for each USB port on your PC – mainly because Windows is not so smart !

2. Install SpectraWiz from appropriate CD folder (Win-9x-ME or Win-2k-NT-XP)

Win2000 only: after installation completes, goto SpectraWiz folder, click on [Fix2000.bat](#).

Note: If folder not found, for USB interface skip to step 3. Otherwise go to website.

3. Install SWUpdate.exe from CD (or website for the most recent SpectraWiz version).

Start SpectraWiz and verify continuous spectral graph display updates appear.

For operation as a SpectroRadiometer, exit SpectraWiz and goto step 4. **If you received a diskette marked "SNS files", copy the files to C:\Program Files\ StellarNet\SpectraWiz.** Otherwise, select the Setup menu to install the wavelength calibration coefficients listed on spectrometer bottom label using "Setup -> Unit Calibration" menu.

4. Install intensity calibration files for SpectroRadiometer operation using the supplied StellarNet diskette. The IRRAD-CAL (irradiance calibration) files are installed by clicking on the file "MyCal-nnnnn.exe", where "nnnnn" = spectrometer serial # shown on label.

Start SpectraWiz and verify continuous spectral graph display updates appear.

Get going FAST -->>> WATCH the SpectraWiz software training videos on CDROM !

Driver and software updates are easily downloaded from the StellarNet website

- | | |
|----------------------------------|---|
| a. All drivers are in SWDrivers | password="wdrivers" |
| b. SpectraWiz for Win95/98/ME | password="wiz" |
| c. SpectraWizNT for WinNT/00/XP | password="wiznt" |
| d. SWUpdate for all Windows vers | password= no password ←get the latest version now |

Additional information is available in the Spectrometer Manual on the StellarNet CDROM.

For PARALLEL cable installation see "Parallel cable installation.doc" on CDROM.

For StellarNet technical support, Phone: 813-855-8687 or email: Support@StellarNet.us