


SPECTRAWIZ UPDATE FOR VISTA USERS

To enable proper use of StellarNet's SpectraWiz software, you must first allow Administrative privileges to the program when using the SpectraWiz software with Windows Vista operating systems.


Instructions:

- 1) Right-click the SpectraWiz program icon
- 2) Select **Properties**
- 3) Click the “**Compatibility**” tab
- 4) Check the “**Run this program as an administrator**” box under the Privilege Level section.
- 5) Click the OK button

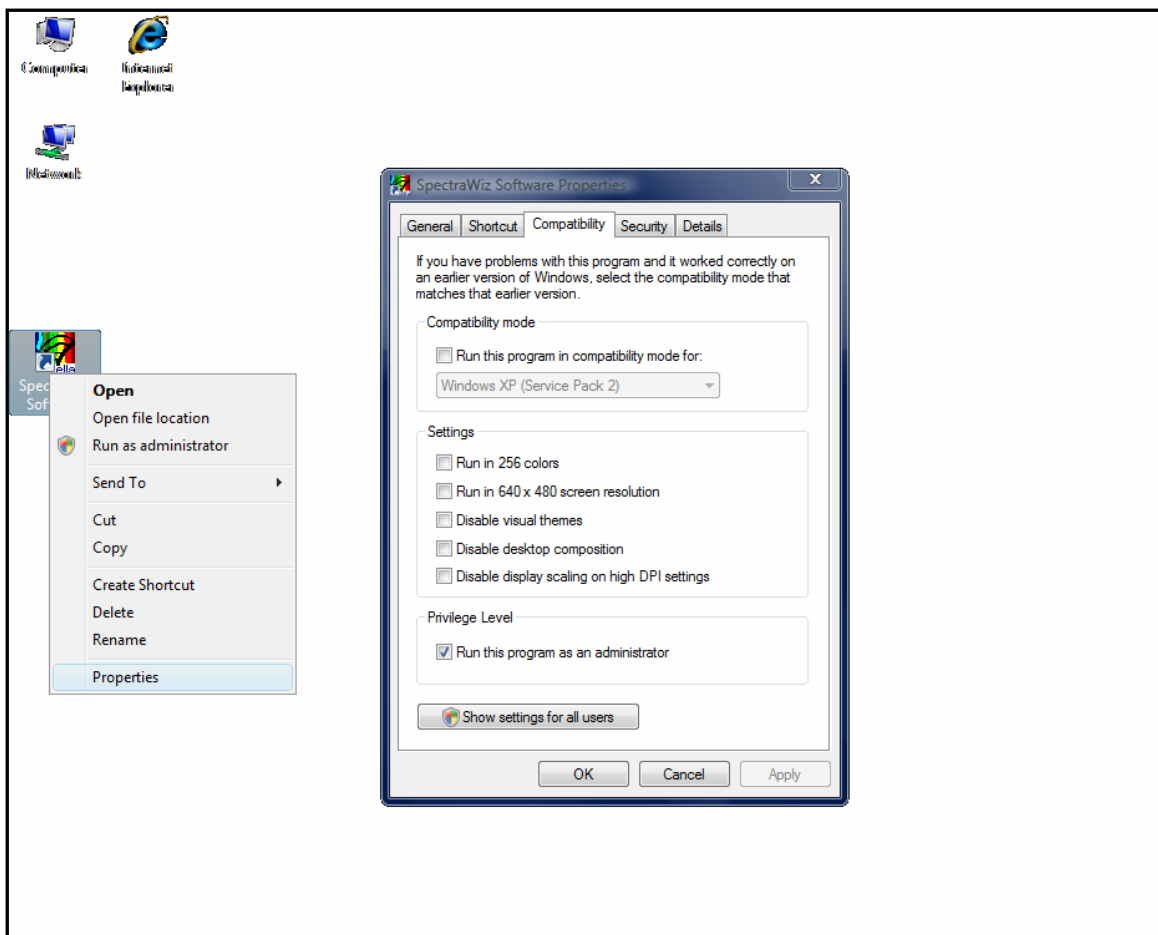


To disable the Windows prompt that asks whether or not to allow access to SpectraWiz every time the software is opened, use the following procedure:

- 1) Double-click the User Accounts icon located in the control panel.

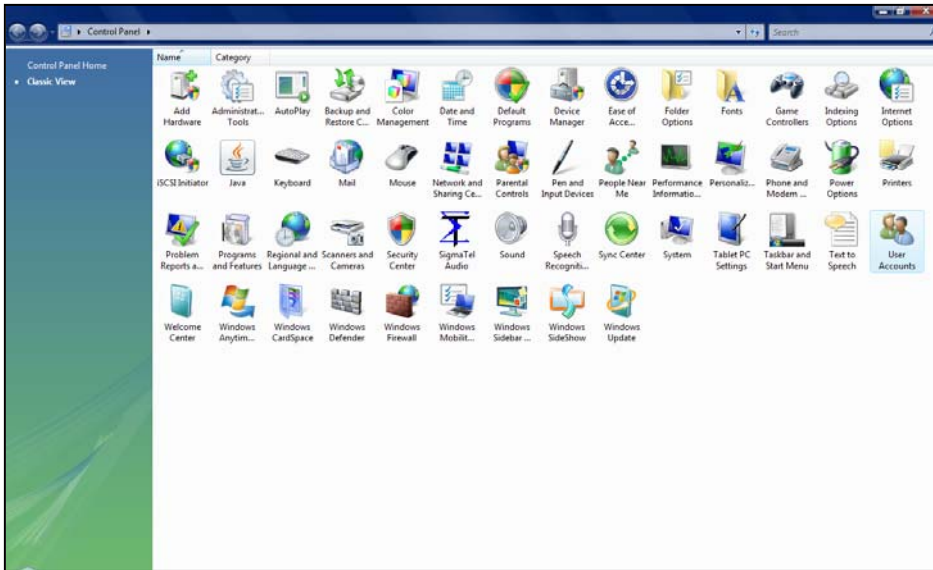

- 2) Select “Turn User Account Control on or off.”

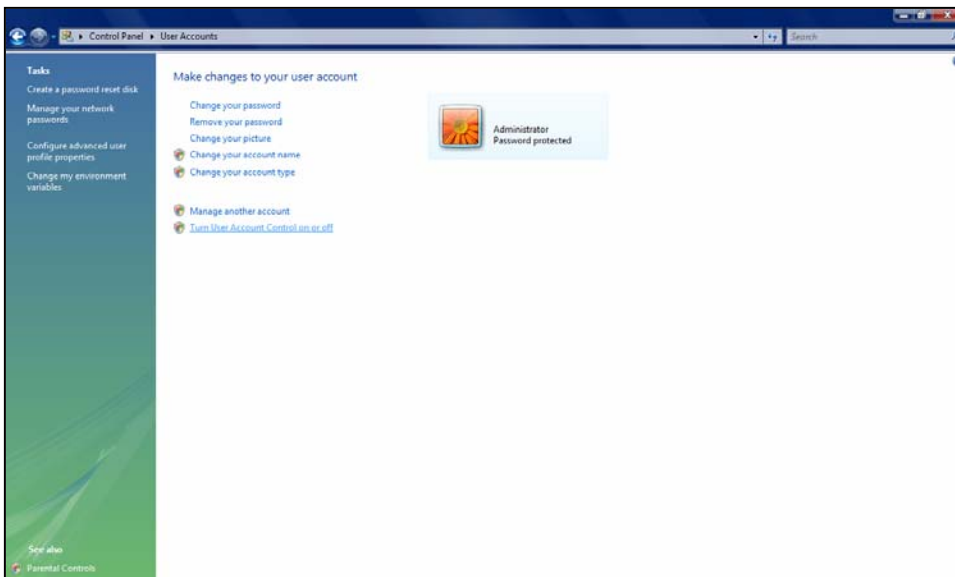

3) Uncheck the box and restart the computer.

